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Undecalcified bone VG stain report**I. Experimental equipment and reagent****Equipment**

Brand name	Manufactures	Model
Oven	Tianjin Labotery instrument equipment Co. Ltd	GFL-230
Microslide	Wanwu	
Optical Microscopy	NIKON	Nikon Eclipse E100
Imaging system	NIKON	Nikon DS-U3

Reagent

Reagent	Manufactures	Cat
Ethanol absolute	Sinopharm Chemical Reagent Co.,Ltd	100092183
Xylenes	Sinopharm Chemical Reagent Co.,Ltd	100023418
VG staining solution	Wanwu	G1046
Neutral balsam	Sinopharm Chemical Reagent Co.,Ltd	10004160
Ethylene glycol ethyl ether acetate	Macklin	E808814-2.5L

II. Experiment procedure**1. Remove the embedding agent and rehydration**

Put sections into ethylene glycol ethyl ether acetate I 6 h at 37 °C and ethylene ether acetate II overnight at 37 °C. Then put them into ethylene glycol ethyl ether acetate III 10-15 minutes at room temperature and ethylene glycol ethyl ether acetateIV 10-15 minutes at room temperature. The sections were rehydrated in 100%I - 100%II - 95% - 90% - 80% alcohol. Each step takes anywhere for 10 minutes. Finally, rinse with running water.

2. Prepare VG staining solution

Mixed with 9 ml of VG staining solution B and 1 mL of VG staining solution A.

3. Washed quickly after use VG staining solution stain for 1 minutes, and three cylinders of ethanol absolute rapidly dehydrate.

4. Then one cylinder of xylenes was used for transparency for 1-5 minutes. Finally, sealed the sections with neutral balsam.

8. Transparent and mount

Microscope observation and collection and analysis images.

III. Result determination

Collagen fibers are red and other tissue components are yellow.

IV. Matter need attention

1. Water should be washed quickly to prevent the red from fading, and ethanol should be removed quickly to prevent the yellow from darkening.