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EDTA antigen retrieval solution (pH 9.0)	Wanwu	G1203
Sodium citrate antigen retrieval solution (pH 6.0)	Wanwu	G1202
PBS solution	Wanwu	G0002
Spontaneous fluorescence quenching reagent	Wanwu	G1221
BSA Primary antibody	Wanwu	G5001
Fluorescent-labelled secondary antibody		
DAPI	Wanwu	G1012
anti-fade mounting medium	Wanwu	G1401

2 Procedure

2.1 Deparaffinize and rehydrate: incubate sections in 2 changes of xylene, 15 min each. Dehydrate in 2 changes of pure ethanol for 5 min, followed by dehydrate in gradient ethanol of 85% and 75% ethanol, respectively, 5 min each. Wash in distilled water.

2.2 Antigen retrieval: immerse the slides in EDTA antigen retrieval buffer (pH 8.0) and maintain at a sub-boiling temperature for 8 min, standing for 8 min and then followed by another sub-boiling temperature for 7 min. Be sure to prevent buffer solution evaporate. Let air cooling. Wash three times with PBS (pH 7.4) in a Rocker device, 5 min each. Use the right antigen retrieval buffer and heat extent according to tissue characteristics.

2.3 Circle and Serum blocking: eliminate obvious liquid, mark the objective tissue with liquid blocker pen. Add 3% BSA to cover the marked tissue to block non-specific binding for 30 min . Cover objective area with 10% donkey serum (for the case of primary antibody originated from goat) or 3% BSA (for the case of primary antibody originated from others)

2.4 Primary antibody: throw away the blocking solution slightly. Incubate slides with primary antibody (diluted with PBS appropriately) overnight at 4 °C, placed in a wet box containing a little water.

2.5 Secondary antibody: wash slides three times with PBS (pH 7.4) in a Rocker device, 5 min each. Then throw away liquid slightly. Cover objective tissue with secondary antibody (appropriately respond to primary antibody in species), incubate at room temperature for 50 min in dark condition.

2.6 DAPI counterstain in nucleus: wash three times with PBS (pH 7.4) in a Rocker device, 5 min each. Then incubate with DAPI solution at room temperature for 10 min, kept in dark place.

2.7 Spontaneous fluorescence quenching : wash three times with PBS (pH 7.4) in a Rocker device, 5 min each . Add spontaneous fluorescence quenching reagent to incubate for 5 min. Wash in running tap water for 10 min.

2.8 Mount: Throw away liquid slightly, then cover slip with anti-fade mounting medium.

2.9 Microscopy detection and collect images by Fluorescent Microscopy. DAPI glows blue by UV excitation wavelength 330-380 nm and emission wavelength 420 nm; FITC glows green by excitation wavelength 465-495 nm and emission wavelength 515-555 nm; CY3 glows red by excitation wavelength 510-560 nm and emission wavelength 590 nm.

3 Results

Nucleus is blue by labeling with DAPI. Positive cells are green or red according to the fluorescent labels used.